
**-ebro-**<sup>®</sup>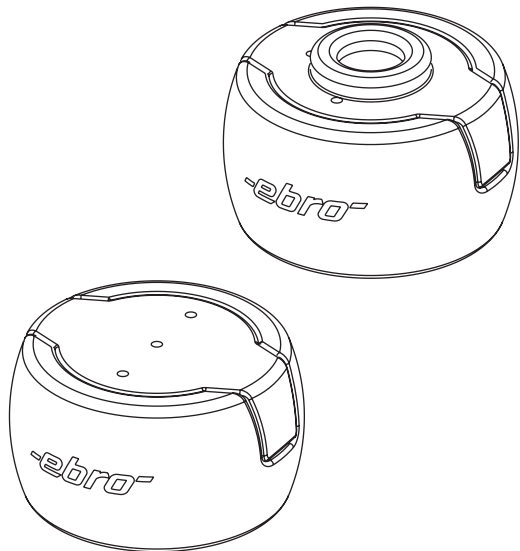

Silikon-Schutzbox  
Silicone protection box  
Boitier de protection en Silicone  
**AL 100/101**

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Please read these operating instructions attentively before operating with your new silicone protection box.

In behave of a safe handling and for protection of the equipment, follow the safety notes marked with the warning sign .

## Safety notes

This box AL 100/101 serves for protection of the data logger series EBI 10 for use in autoclaves and/or for use at higher temperatures.

This box is not a thermal protection box in the conventional sense.

**Take care when using the box at high temperatures! Note when taking the Silicone protection box with the logger from the autoclave it is very hot.**

**Do not grasp the logger with bare hands immediately after use! Danger of burning !**

**Please remove the data logger from the thermal insulation box immediately after the measurement is completed, in order to prevent possible heat damage of the logger due to the thermal inertia of the insulation box.**

Follow the diagram and the technical data on the next page also the safety notes described before to this!

**Never exceed the maximum permitted temperature of 150°C!**

**Take care when opening the protection box as the logger can be very hot.**

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